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Filterwechsel UFA.20.30+ / UFA.20.30*

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1. NOTES - GUIDELINES

1.1. SAFETY NOTES

- Take notice of the legislation and provisions to be applied in dealing with hazardous materials.
- Only use the recirculating air filter system in a proper condition.
- Observe all notes in these instructions.
- Follow the instructions of the safety specialists, technical monitoring services (BG, GAA, pharmaceuticals councils), construction supervision etc.
- Observe accident prevention regulations and workplaces ordinances.
- The filter cannot be washed out. It must always be replaced.
- Protect yourself when changing the filter by wearing personal protective equipment (e.g. safety goggles, respiratory protection, safety gloves etc.).
- Change the filter in a safe place (e.g. hazardous material workplace etc.).

1.2. GENERAL NOTES

These operating instructions are for practical use and should be available to the user where the recirculating air filter system is to be used. Keep these operating instructions best of all next to the filter attachment. An assured, faultless functioning of the recirculating air filter system is only ensured when these instructions are followed.



CAUTION:

The recirculating air filter system is only to be repaired by specialists trained for this very purpose. Given damage the appliance is to be repaired or replaced by the manufacturer. Since the filter is contaminated with pollutants, it is to be replaced only by the manufacturer's trained personnel or authorised persons.

After replacement, the contaminated active carbon filter must be properly disposed of or submitted for recycling.

Development and manufacture: asecos GmbH, Weiherfeldsiedlung 16-18, D-63584 Gründau
Serial number: Please take the serial number from the type plate

2. CHANGING THE FILTER IN THE UFA.20.30+

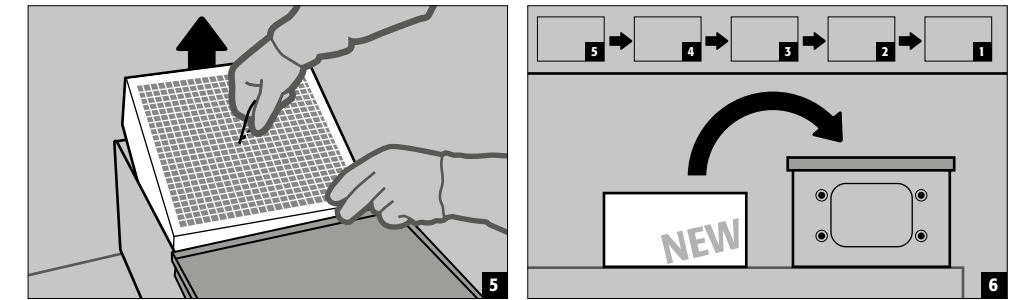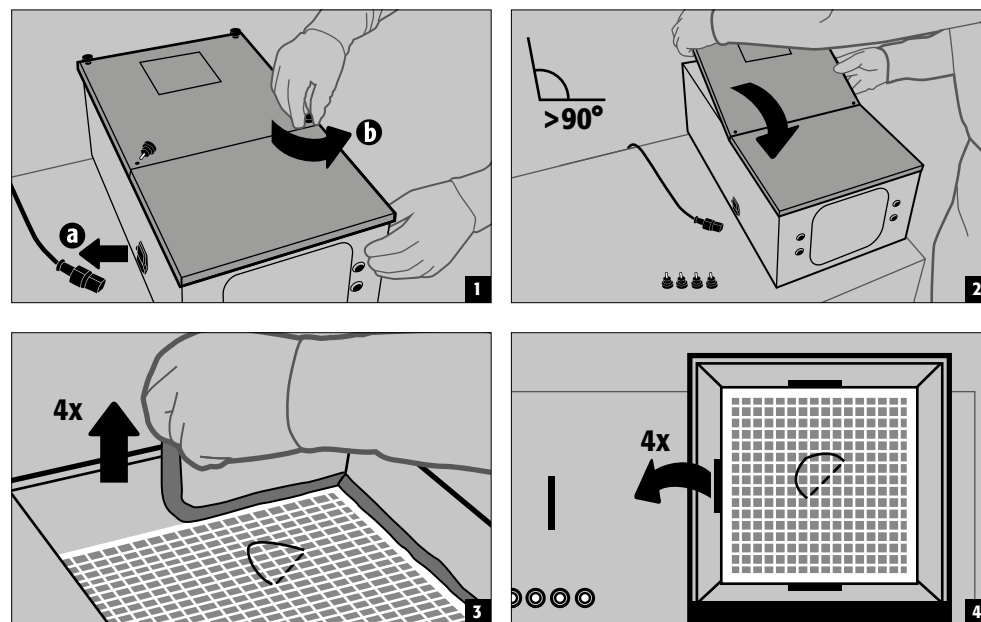CAUTION:

The filter must be disposed of properly or taken to an official recycling collection point. Returned filters will be disposed of by the manufacturer. Once the filter, centring pieces and sealing cords have been fitted, RESET the filter saturation monitor. ► SEE 3

3. RESET THE FILTER SATURATION MONITOR IN THE UFA.20.30+


CAUTION:

The filter saturation monitor must be reset after each filter change

SERVICE: In case of questions please contact Customer Service 01805 92 20 92* (Germany) or +49 6051 9220-0 (foreign countries).

* 14 ct/ per minute or part of a minute from the landline network of the Deutsche Telekom AG.
Max. 42 cent/minute from mobile phones.

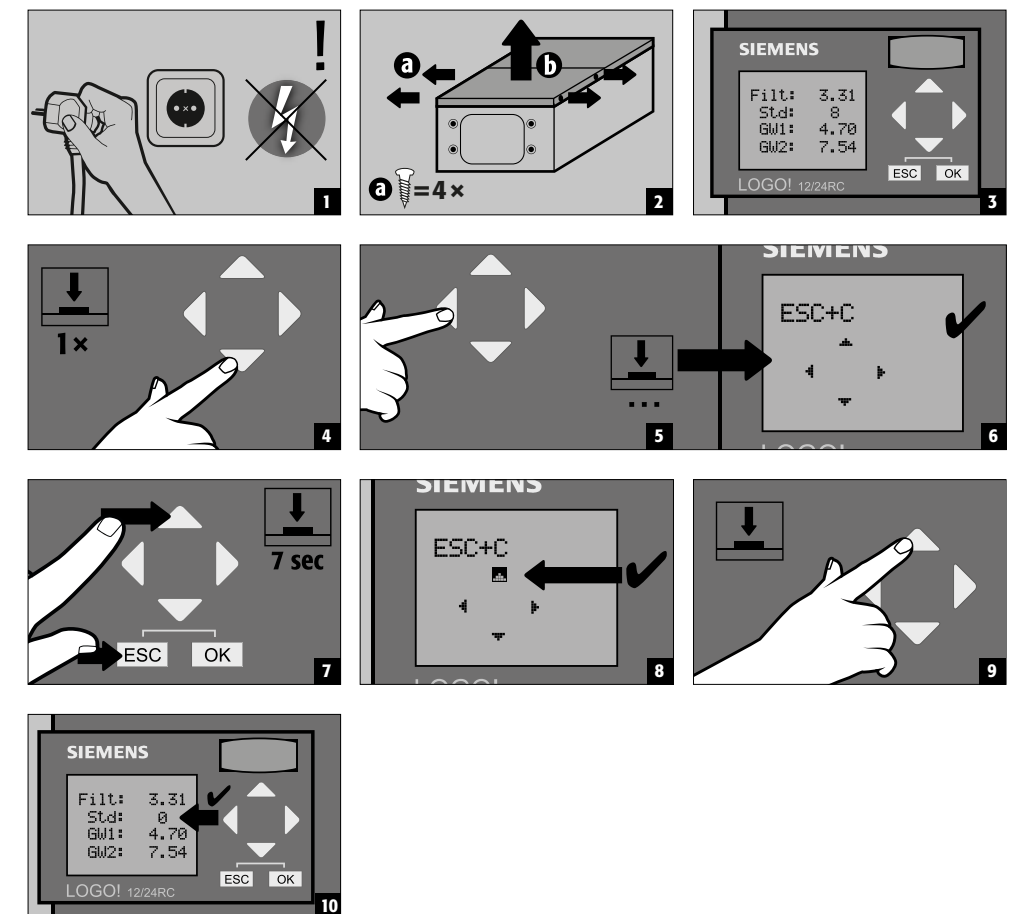4. CHANGING THE FILTER IN THE UFA.20.30*










CAUTION: General safety notes and guarantee

The UFA.20.30* recirculation filter is a safety-relevant equipment and must be checked at least once per year, or more frequently depending on the level of use. Inspections and filter changes must be carried out by specialised staff authorised by the manufacturer, using original spare parts from the manufacturer.

If this inspection is not carried, or not carried out by specialised staff authorised by the manufacturer, the customer's guarantee claims against the manufacturer will be invalidated.

Protective measures

If the recirculating air filter system is used as intended, all air conducting parts will be contaminated. Therefore additional protective measures such as the use of personal protective equipment are required when changing the filter.

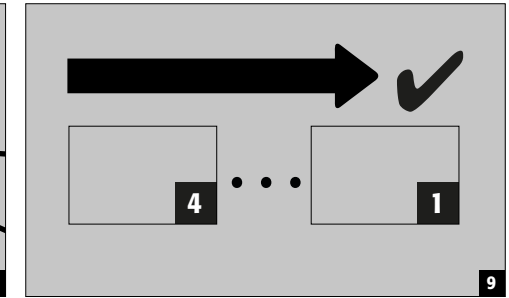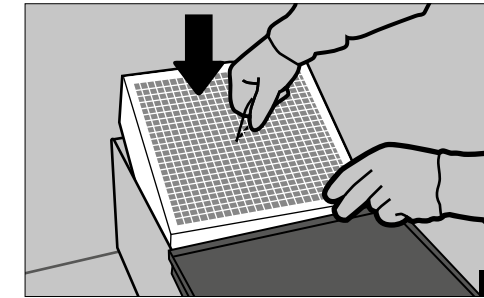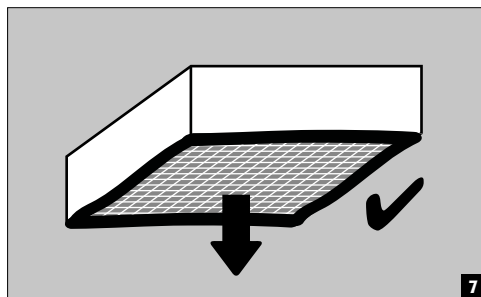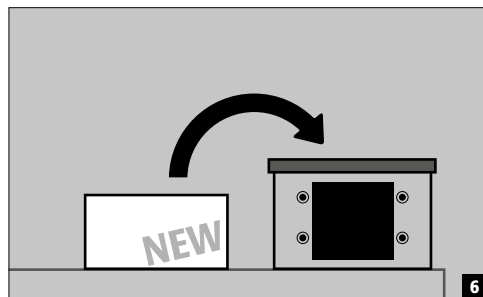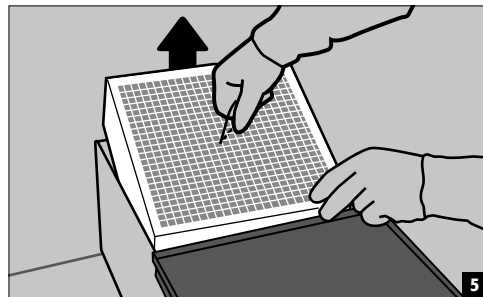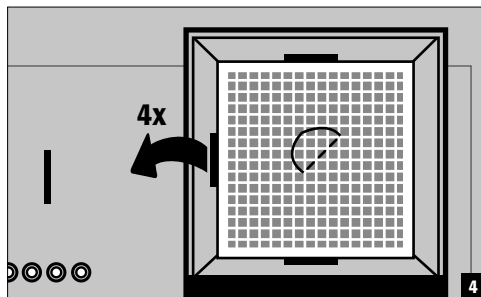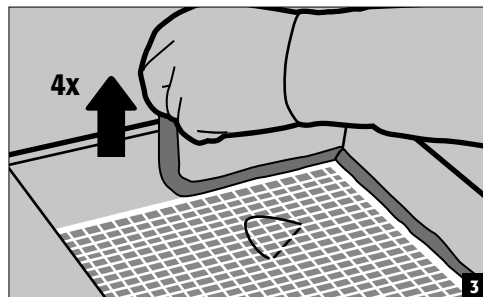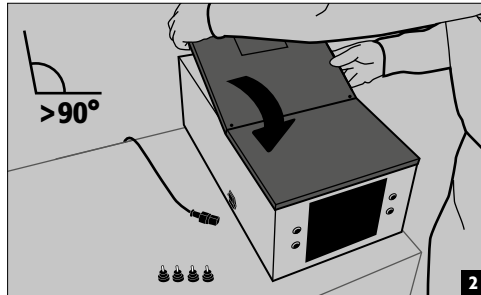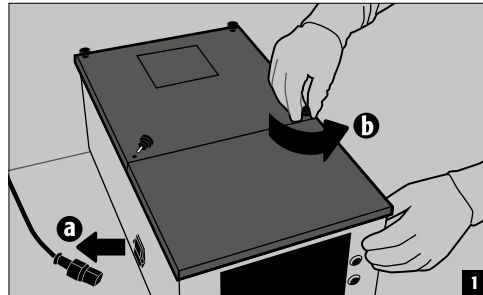CAUTION:

Make sure that the gasket does not touch the housing (figure 7), in particular that it does not touch the rear upper bar. In order to prevent damage to the cables, the filter must not touch them. Check whether cables or lines are damaged. Check that the protective grounding wire is correctly connected to the cover.

5. RESET THE FILTER SATURATION MONITOR IN THE UFA.20.30*

CAUTION:













The filter saturation monitor must be reset after each filter change.

SERVICE: In case of questions please contact Customer Service 01805 92 20 92* (Germany) or +49 6051 9220-0 (foreign countries).


* 14 ct/ per minute or part of a minute from the landline network of the Deutsche Telekom AG. Max. 42 cent/minute from mobile phones.

5.1. MODELS BEFORE 2014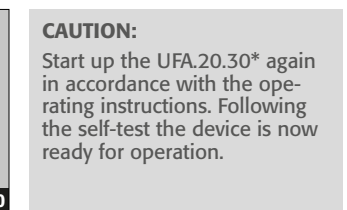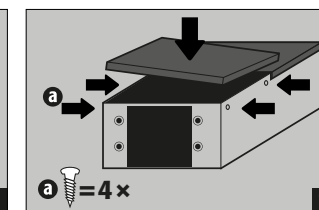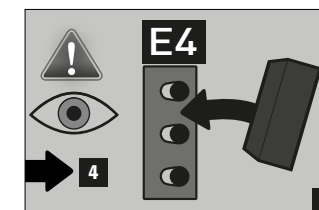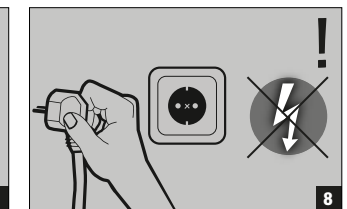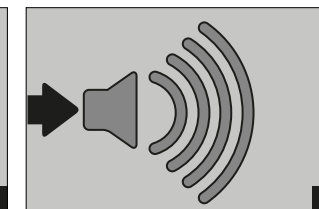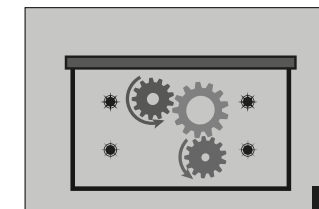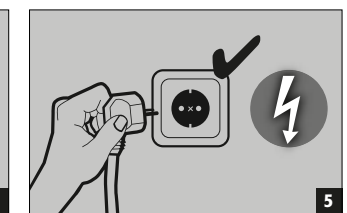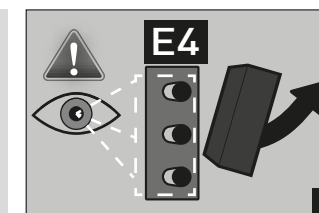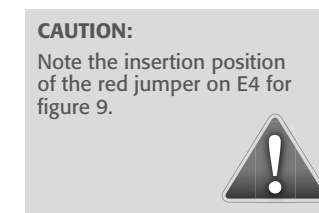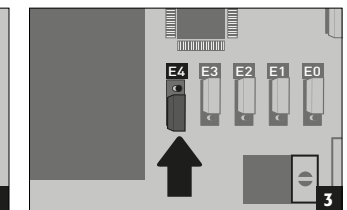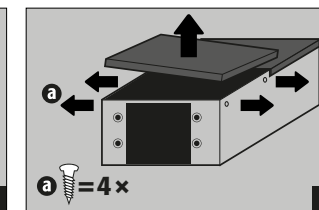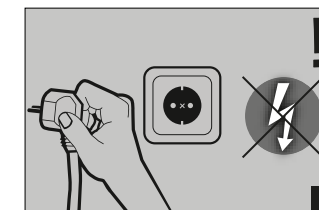


5.2. MODELS FROM 2014

Resetting of the circuitry by persons with no knowledge of electronics: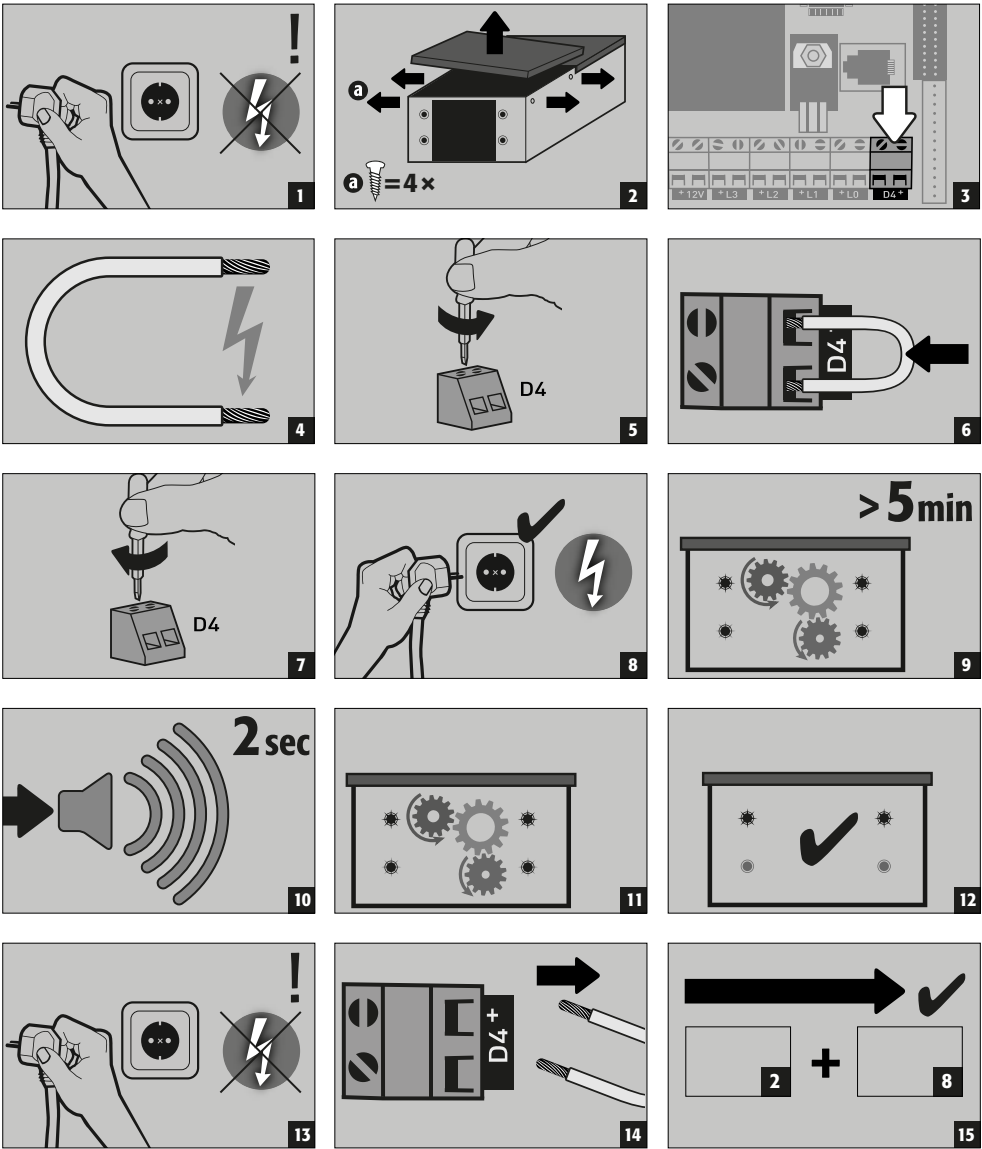


CAUTION
Start up the UFA.20.30* again in accordance with the operating instructions. Following the self-test the device is now ready for operation.

Resetting of the circuitry by an electrical specialist: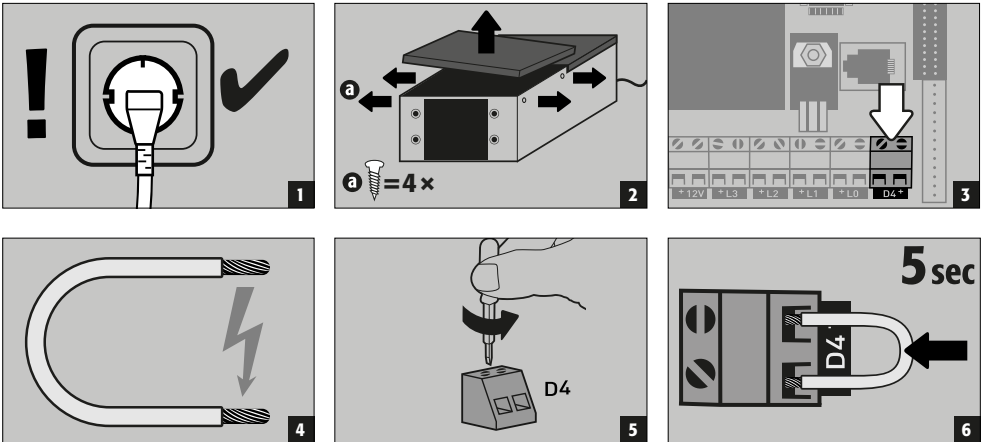

9 self test

11 warm-up period

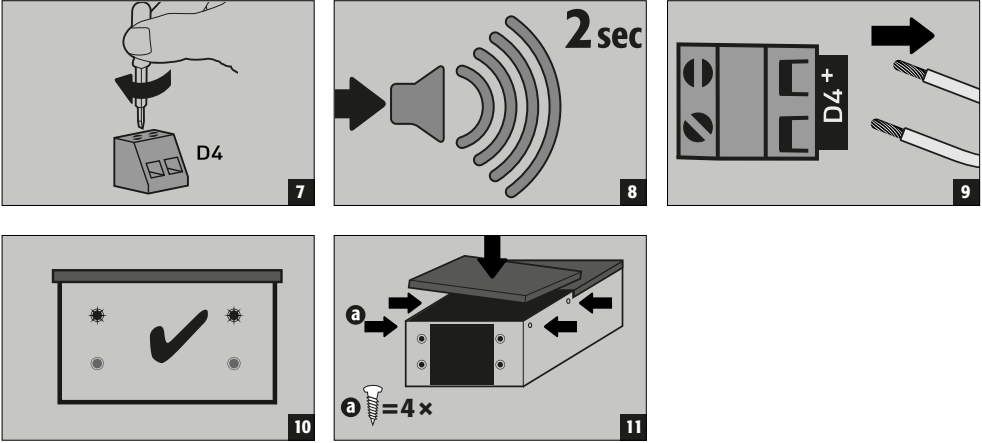CAUTION:
Following the self-test the device is now ready for operation.

6. FUNCTIONAL CHECK AND CARE

To ensure reliability, the recirculating air filter system should be serviced and checked as to functioning by a specialist at least once a year. At the same time the fitted filter unit is to be replaced at least once a year - or sooner should the saturation limit be reached. The integrated ventilation and filter saturation monitoring system electronically checks on functionality. As a result, the recirculating air filter system function is constantly examined during operation.

The activated carbon filter of the recirculating air filter system must be replaced soon after the red LED of the filter display blinks. The filter can be changed by one of our trained service technicians and must be requested promptly by calling the service hotline on 0180 - 5 92 20 92* (Germany) or +49 6051 9220-0 (foreign countries).

Having the filter changed by asecos service personnel ensures that contaminated activated carbon is recycled in a defined disposal chain. Clean the outside of the recirculating air filter system with a soft cloth and a common, mild household cleaner.

CAUTION:
Make sure that no moisture can penetrate into the inside of the recirculating air filter system during cleaning.

SERVICE: In case of questions please contact Customer Service 01805 92 20 92* (Germany) or +49 6051 9220-0 (foreign countries).
* 14 ct/ per minute or part of a minute from the landline network of the Deutsche Telekom AG. Max. 42 cent/minute from mobile phones.

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* 14 centesimi / per ogni minuto iniziato dalla rete fissa della Deutsche Telekom AG.