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Sicherheitsschränke V-LINE

## USER MANUAL

Safety storage cabinets V-LINE

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### 10 YEARS WARRANTY\*

\* Upon conclusion of an asecos service and maintenance agreement (**PREMIUM tariff**), you will get a warranty extension for up to 10 years for your V-LINE safety storage cabinet.

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For detailed information about our warranties please visit [www.asecos.com/service-conditions](http://www.asecos.com/service-conditions)

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asecos GmbH

## **YOUR PERSONAL DOCUMENTATION TO THE asecos SAFETY CABINET**

Dear Customer,

you have made a decisive investment in safety for your company by purchasing this asecos safety storage cabinet. You now own an innovative product made of top-quality materials guaranteeing the highest quality standards.

asecos safety storage cabinets have complete authorisation documents. We archive the authorisation documents for every individual cabinet, keeping them ready for you should you ever need them (e.g. for a works inspection or similar). Simply request them using this form.

Tear off or copy that page and return to us by fax with your address and serial number of the cabinet on it.

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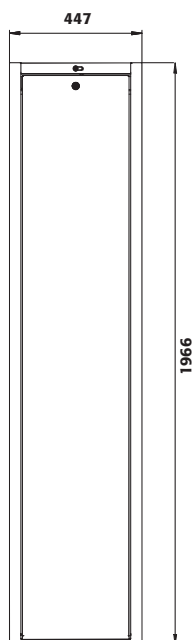
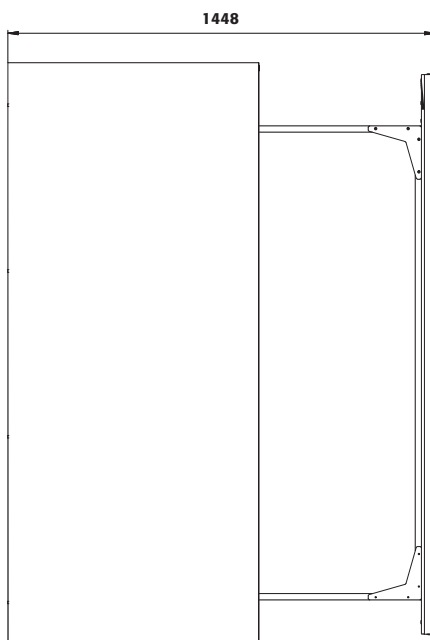
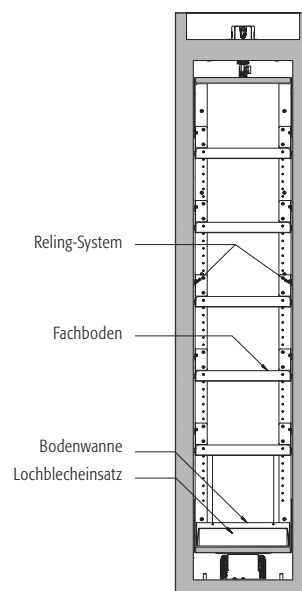
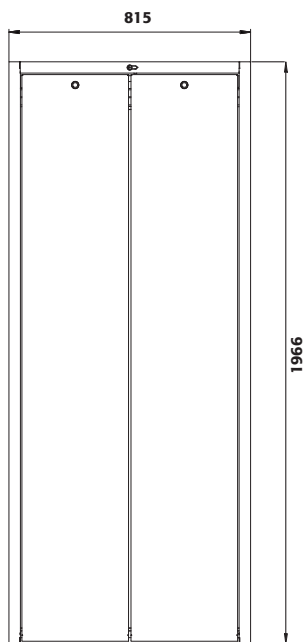
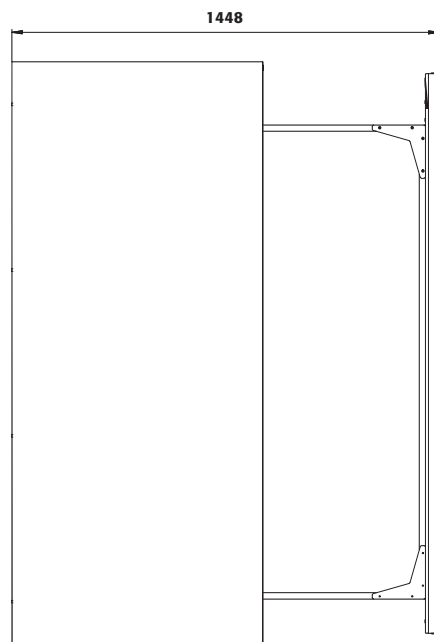
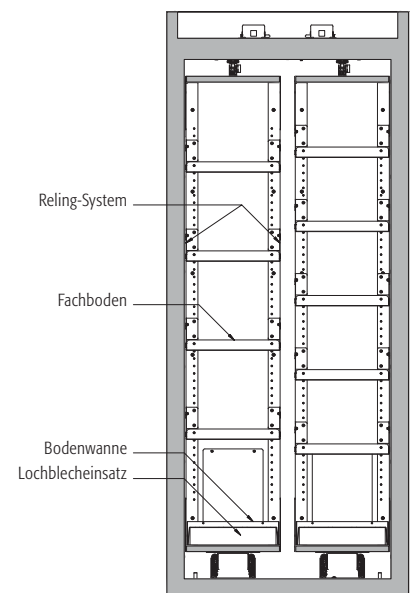
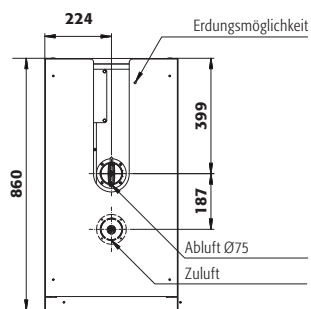
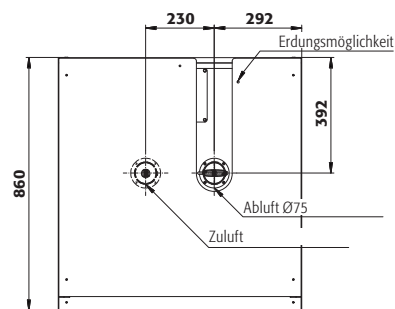
# OPERATING INSTRUCTIONS

Dear customer,

Thank you very much for purchasing a safety storage cabinet from our company, with which you have made a decisive investment in the safety within your company. Our safety storage cabinets make the storage of hazardous materials at the workplace safe and convenient for you.

Please read these operating instructions very carefully. Get to know the advantages and simple operability of our safety storage cabinets in detail. This simplifies the daily handling of hazardous materials for you.

Many thanks.  
Your asecos team

---

## CONTENT

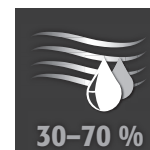
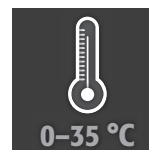
<b>1. NOTES ▪ GUIDELINES ▪ GUARANTEE</b>	<b>20</b>
1.1. general safety notes	20
1.2. guarantee	20
1.3. Cabinet Details	20
<b>2. TRANSPORT</b>	<b>21</b>
2.1. Dismantling of the transport packaging	21
2.2. In-plant transport	22
<b>3. ERECTION ▪ COMMISSIONING ▪ FUNCTION</b>	<b>22</b>
3.1. commissioning	22
3.2. Alignment Of The Cabinets	22
<b>4. CLOSING MECHANISMS</b>	<b>23</b>
4.1. locking of the cabinet	23
4.2. Unlocking of the cabinet	23
4.3. Automatic door closing system (TSA)	23
4.4. disable the automatic door closing system	24
<b>5. SAFETY FUNCTIONS</b>	<b>24</b>
5.1. Closing the drawer during blackout	24
5.2. Automatic closing in the event of fire	25
5.3. Blocking the drawer while opening	25
5.4. Blocking the drawer while closing	25
<b>6. INTERIOR FITTINGS</b>	<b>26</b>
6.1. Spill Containment Sump	26
6.2. Shelves (Height-Adjustable)	26
<b>7. STORAGE</b>	<b>27</b>
<b>8. EARTHING</b>	<b>27</b>
<b>9. VENTILATION</b>	<b>27</b>
<b>10. FIRE ▪ DISPOSAL</b>	<b>28</b>
10.1. Opening The Cabinet After The Fire	28
10.2. Disposal	28
<b>11. SAFETY CHECKS</b>	<b>28</b>
11.1. In general	28
11.2. Battery check	29
<b>1. TECHNICAL DATA</b>	<b>30</b>

## 1. NOTES - GUIDELINES - GUARANTEE

### 1.1. GENERAL SAFETY NOTES

- Observe applicable statutes and regulations, and the notes in these operating instructions, when handling hazardous materials
- Work on the electrical system is to be carried out only with the power turned off and only by qualified electricians – refer here to the regulations of the local electricity supply company.
- The on-site installation conditions are to be observed (e.g. bolting the cabinets to the building).
- The instructions of the supervisory engineering department must be followed.
- Observe accident prevention regulations and workplace ordinance
- Ensure that the necessary safety checks are only carried out by authorised staff using original spare parts
- Only use the cabinet after having been properly instructed; access is to be forbidden to unauthorised persons.
- The pivoting area of the doors is to be kept free at all times; doors/drawers are to be kept closed
- By assigning trained/authorised technical personnel you can prevent the malfunctions, damage and corrosion damage that result from inappropriate transport.
- Observe the upper limits for stored quantities, loading etc.
- Containers with a volume larger than that of the bottom collecting sump may not be placed inside the cabinet; spilt hazardous materials are to be collected immediately and removed
- Vessels containing aggressive chemicals (acids/alkalis) are to be kept in special cabinets for acids and alkalis or on shelves for acids and alkalis as well as in cabinets with metal-free interior equipment.
- Before storing such materials, check that the cabinet's surface is resistant to the chemicals.
- Before the initial commissioning, the safety storage cabinet is to be examined by the user for possible damage.

#### Set-up requirements

### 1.2. GUARANTEE

The guarantee for this product is agreed between you (the customer) and your dealer (the seller). As the manufacturer, asecos guarantees the products listed in the operating instructions for a period of 24 months from the date of delivery. All model safety equipment are subject to a compulsory annual inspection by specialised staff authorised by the manufacturer. Otherwise the customer's guarantee claim against the manufacturer expires.

### 1.3. CABINET DETAILS

A complete overview of the models can be found at the start of the operating instructions.

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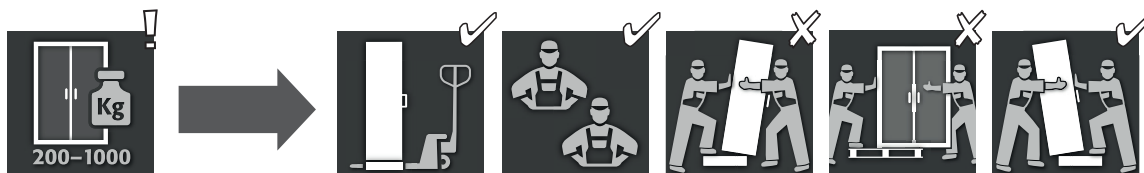
Cabinet data Logbook (included with the cabinet)  
Technical drawing Appendix 1  
Technical data Table in Appendix 2

#### V-LINE

These models are tested, certified and marked according to requirements of BS EN 14470-1. They are intended for the storage of flammable liquids in work rooms in accordance with the valid national regulations.

## 2. TRANSPORT

**CAUTION:** Transport the cabinet in an upright position on a pallet truck, tied and secured against slipping, until the final place of installation is reached. The transport locks in the door joints may only be removed directly at the place of installation! Inappropriate transport can lead to concealed damage to the fire protection insulation! We can only guarantee the necessary quality if the cabinet is transported to the place of its use by our specially trained staff.


**CAUTION:**  
The cabinet must not be moved by pulling the opened drawer! Pulling or sliding may only be carried out in the body housing!

### Tilting the cabinet (may only be done without jolts)
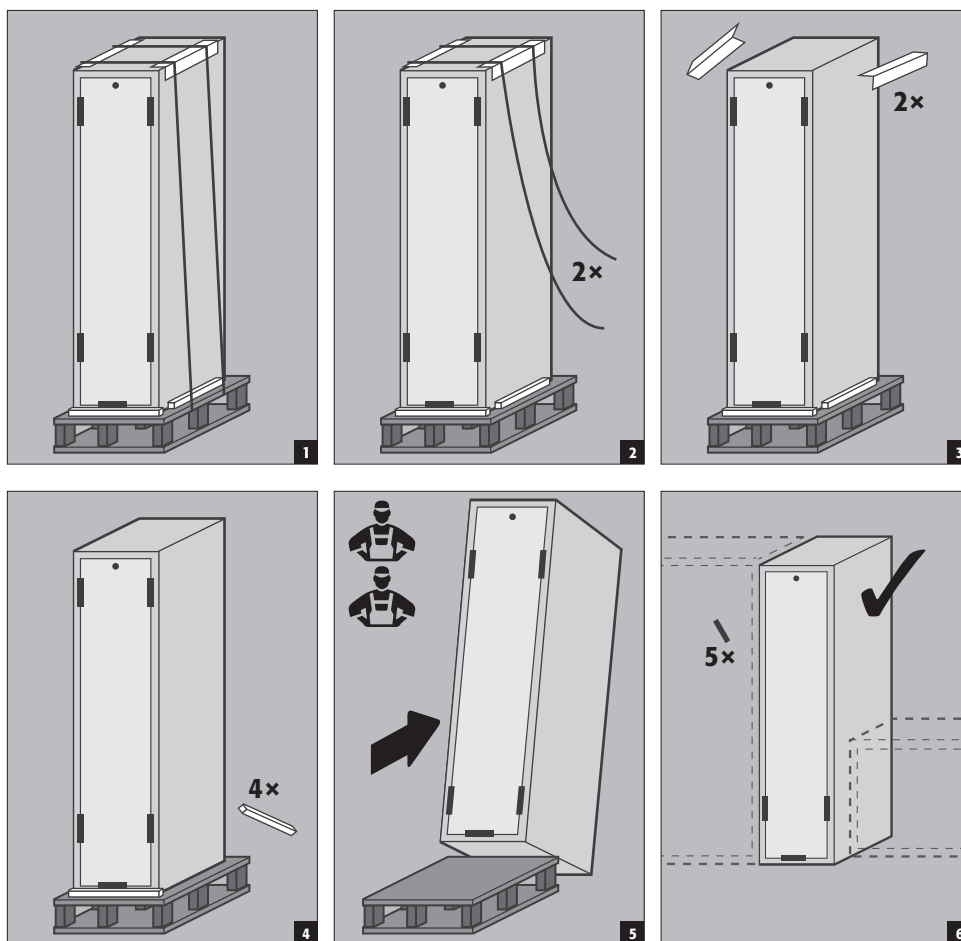
### 2.1. DISMANTLING OF THE TRANSPORT PACKAGING


## 2.2. IN-PLANT TRANSPORT

### CAUTION:

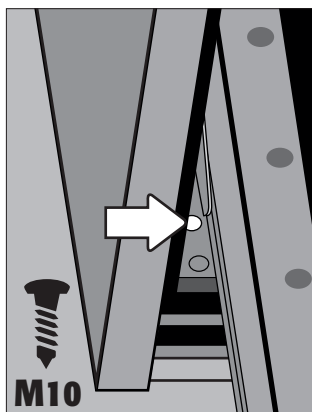The cabinet must not be moved by pulling the opened drawer! Pulling or sliding may only be carried out in the body housing!

### Steps for manoeuvring the cabinet:

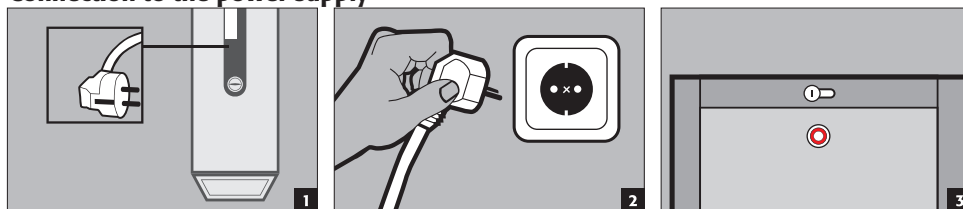
- Open the drawer of the cabinet
- Connect the belts with steel angles or steel rings using a screw M10
- Screw the steel angles/rings into the threaded insert in the base of the cabinet (pics 1 and 2)
- Move the cabinet carefully using the belts

## 3. ERECTION - COMMISSIONING - FUNCTION

### 3.1. COMMISSIONING

Before putting into operation for the first time, the user must carry out an examination of the safety storage cabinet for possible damage, such as defective or loose sealing elements, correct alignment and perfect functioning of the door elements. Use the cabinet and accessories only if they are in an orderly condition.

#### Connection to the power supply


### CAUTION


The cabinet must be connected to the power supply at least 1 hour before filling it the first time!

### 3.2. ALIGNMENT OF THE CABINETS

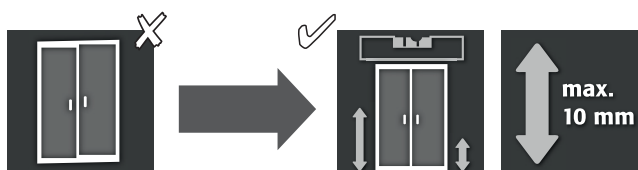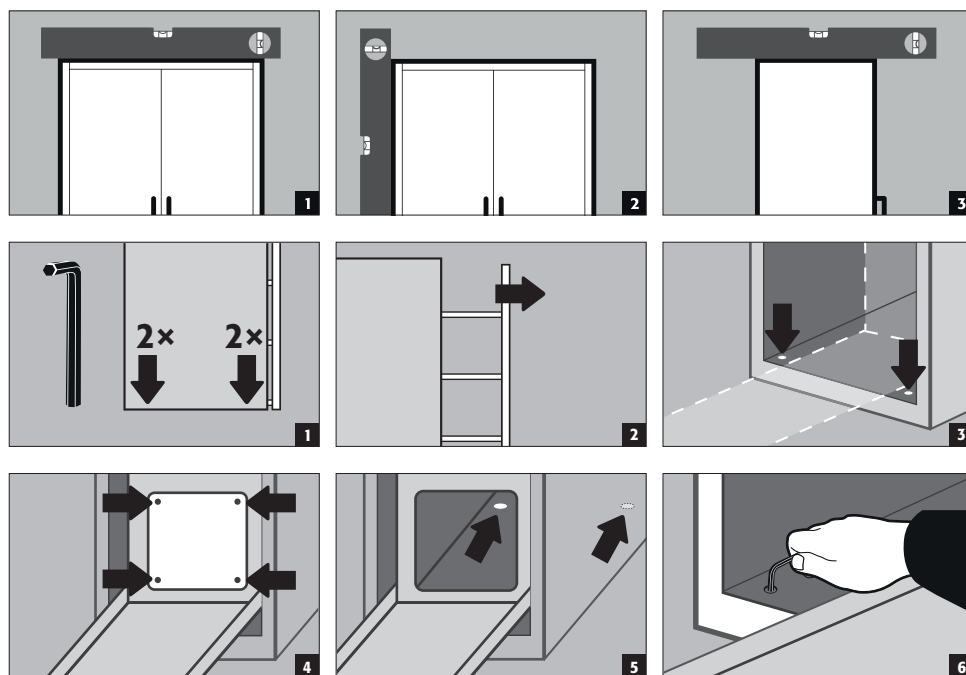
After disabling the automatic door closing (see 4.4) and opening the drawer, the cabinet can be aligned using the integrated adjusting aids.


## 4. CLOSING MECHANISMS

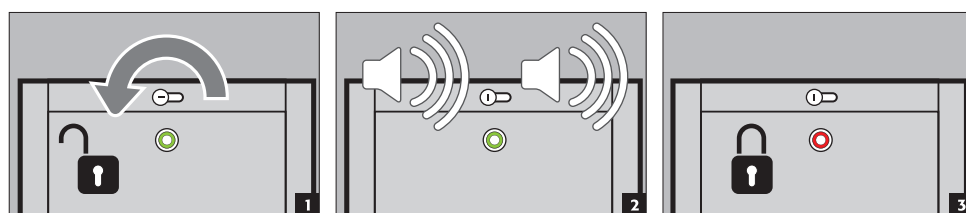
### 4.1. LOCKING OF THE CABINET


### 4.2. UNLOCKING OF THE CABINET

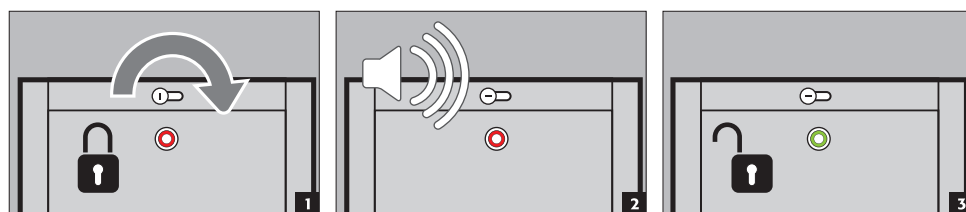
### 4.3. AUTOMATIC DOOR CLOSING SYSTEM (TSA)
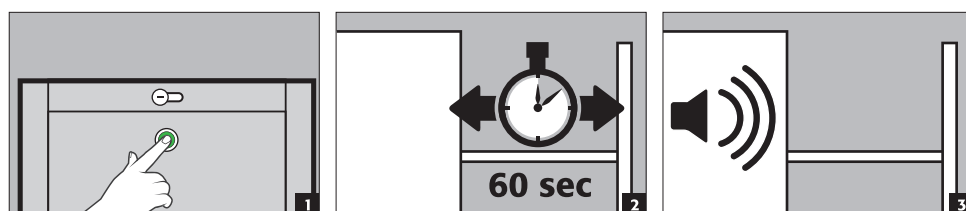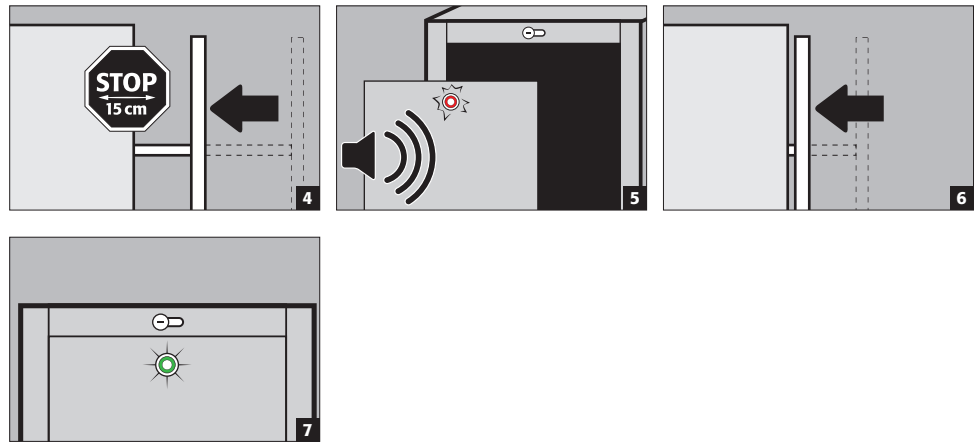
#### 4.4. DISABLE THE AUTOMATIC DOOR CLOSING SYSTEM
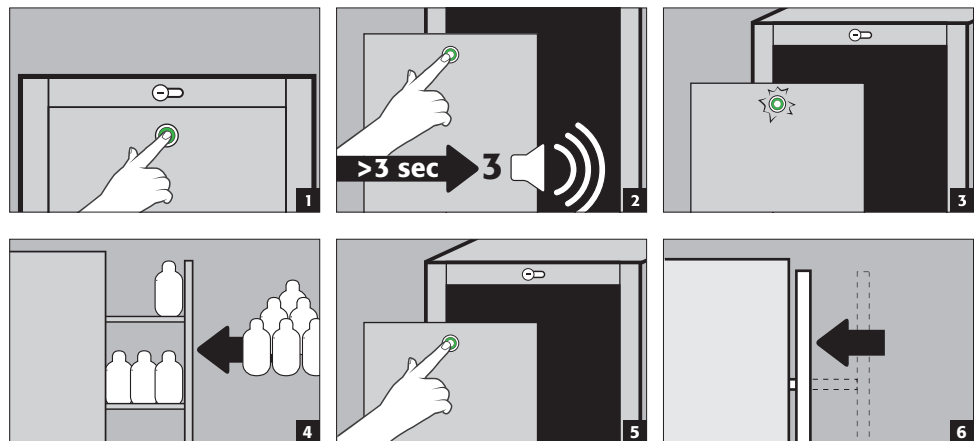
#### ATTENTION:

In case of fire the automatic door closing is always ensured!

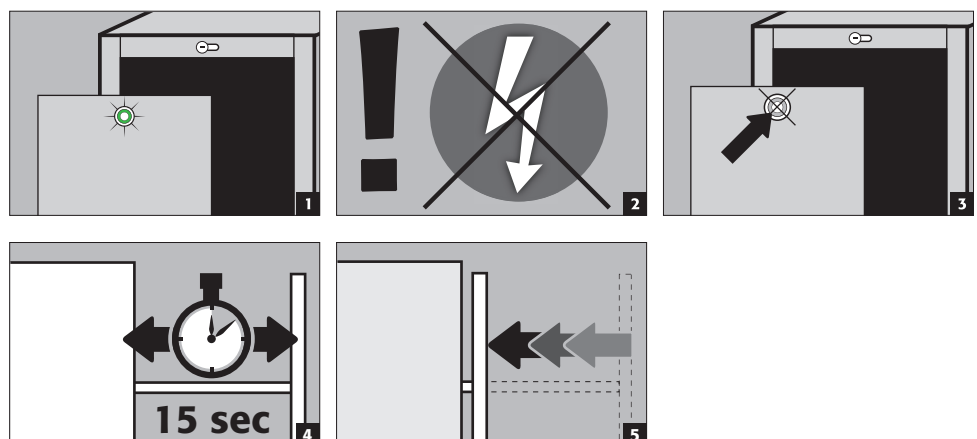
## 5. SAFETY FUNCTIONS

### 5.1. CLOSING THE DRAWER DURING BLACKOUT


#### CAUTION:


During a blackout the opened drawer will close automatically using an integrated battery. Please note, it only has power **for one** closing routine.

## 5.2. AUTOMATIC CLOSING IN THE EVENT OF FIRE

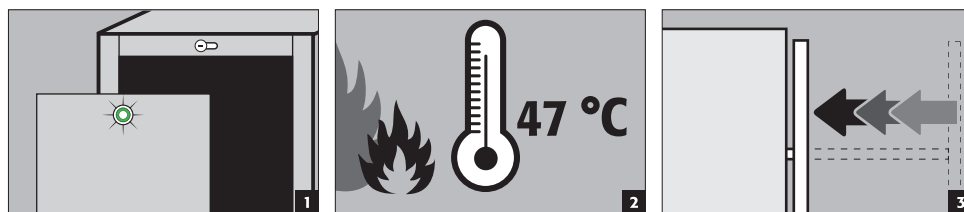
## 5.3. BLOCKING THE DRAWER WHILE OPENING
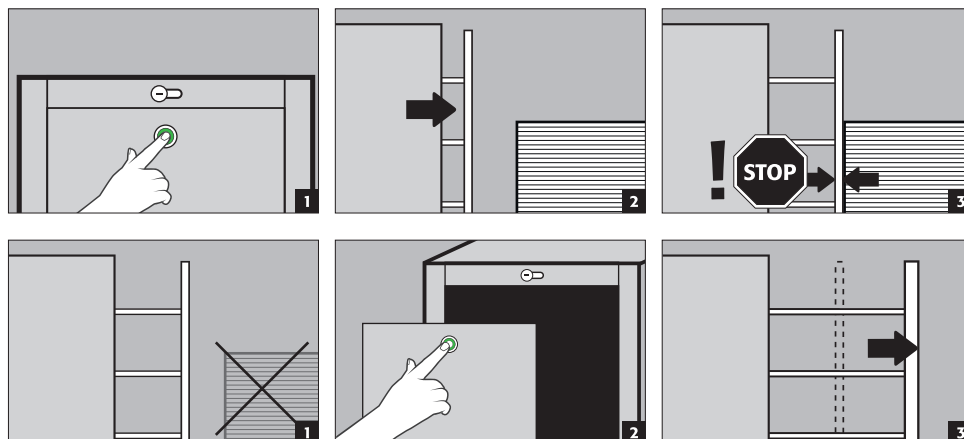
### CAUTION

If no direct user intervention occurs after the blockage, the cabinet will give an acoustic and visual alarm after 20 seconds. In that case, please remove the blockade and follow the previous steps!
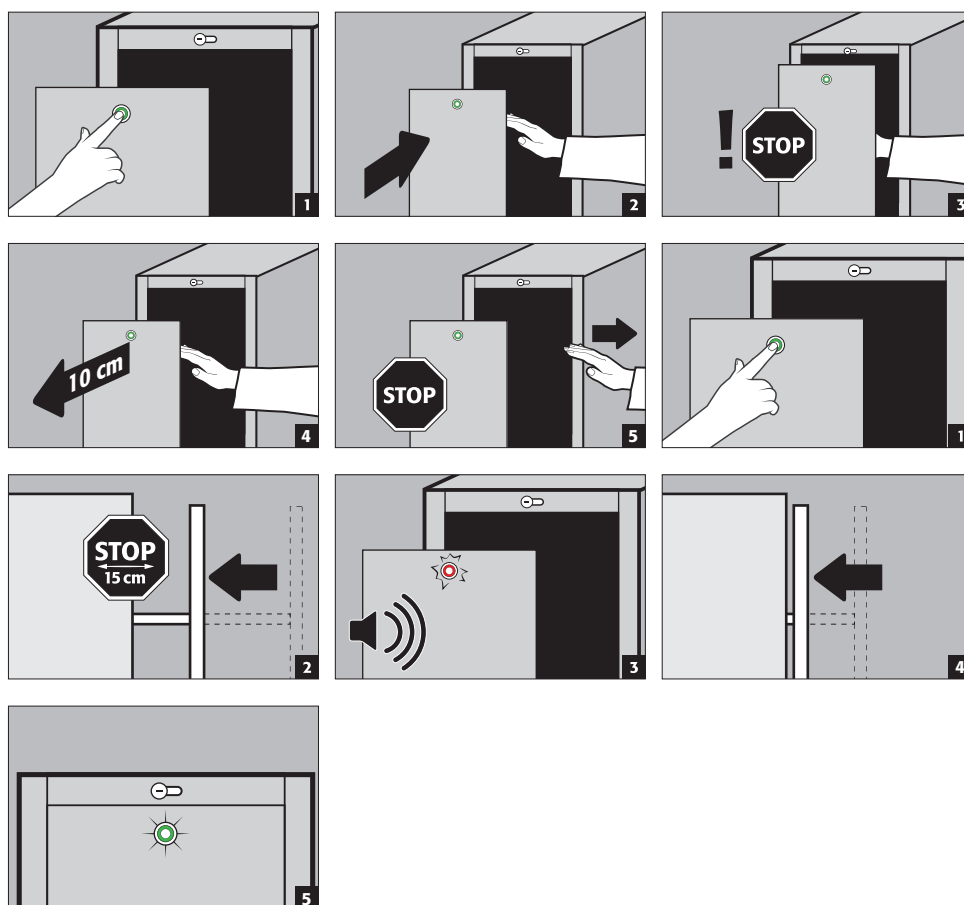
## 5.4. BLOCKING THE DRAWER WHILE COSING

### CAUTION

If no direct user intervention occurs after the blockage, the cabinet will give an acoustic and visual alarm after 20 seconds. In that case, please remove the blockade and follow the previous steps!

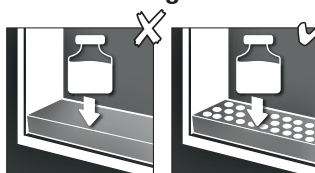
## 6. INTERIOR FITTINGS

### 6.1. SPILL CONTAINMENT SUMP

According to EN 14470-1: A spill containment sump must be installed below the lowest storage level. The spill containment sump must be able to hold a minimum volume of 10 % of all the containers stored in the cabinet [1] or at least 110 % of the volume of the largest single container [2] , whichever of these volumes is larger.

**Use as a storage surface in case of cabinets with several storage levels:**



### CAUTION:




Use as a storage surface is possible only in conjunction with a perforated metal plate insert

### Leaks:

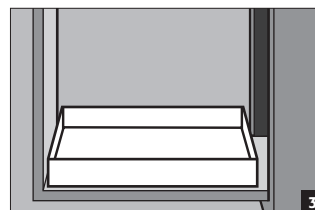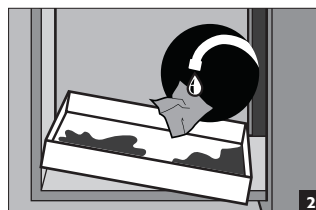- Liquid in the sump is to be collected using suitable means. The choice of means is your own responsibility.

### 6.2. SHELVES (HEIGHT-ADJUSTABLE)
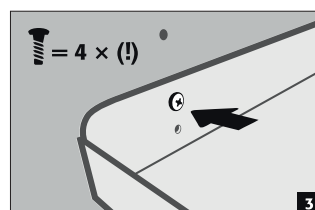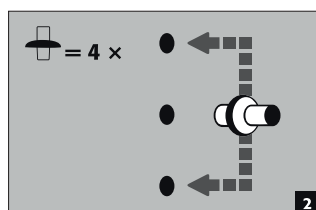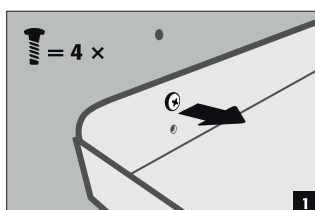
## 7. STORAGE

- Vessels containing aggressive chemicals (acids and alkalis) are to be kept in special cabinets for acids and alkalis or on shelves for acids and alkalis as well as in safety storage cabinets with metal-free interior equipment.
- The storage of corrosive liquids can affect the function of the shut-off equipment for the air supply and exhaust.
- Each container placed inside reduces the minimum containment volume in relation to the total stored quantity.

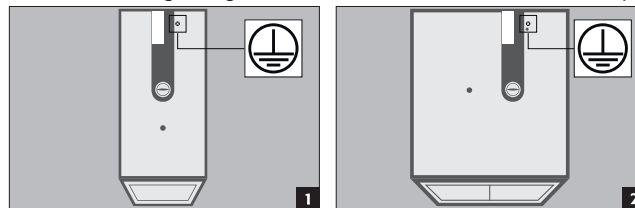
Applicable for all models:  
Depending on the type class of the cabinet, observe the respective valid national regulations!

## 8. EARTHING

An earth connection may be necessary, depending on the customer-specific use of the safety storage cabinet.

The decision regarding this is taken on the basis of a risk analysis by the customer.

### Earthing when decanting:

Connection to a potential equalisation system is compulsory in this case. Cabinets in which decanting can be performed should additionally be equipped with an earth terminal in the interior (at a surcharge on enquiry) which must be connected to the decanting containers. This is standard with cabinets from the UB-LINE with drawers.

The spare parts required for the connection are not part of the scope of supply.

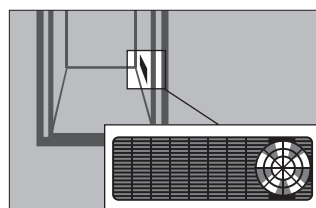
## 9. VENTILATION

**CAUTION:** The fire protection valves near the supply and extraction air connections are safety and maintenance-relevant components. In order to check the correct operation or to replace a defective element, please ensure that flexible connecting hoses or sliding pieces are used to connect to an extraction system in order to ensure simple dismantling and re-mounting of the extraction air connection.

- All models can be connected to a technical exhaust system via the exhaust connector. Please see the respectively applicable national regulations for the specified air recirculation rate, maximum pressure losses and exhaust air ducts.
- If installed on an exhaust system, a check must be carried out to ensure proper connection to the system (e.g. using a smoke tube)

### Ventilation wheel for technical ventilation

- The ventilation wheel serves as an indicator for correct ventilation. ▪ If the wheel turns, the necessary 10-fold air recirculation rate is exceeded.

**CAUTION:** In general these models can be operated without technical ventilation. The direct environment of the cabinet can be an explosive zone in this case. If the cabinet is to be operated without an exhaust connection, the user must attach a marking to it. With some models the supply and exhaust connectors are not pre-mounted. They are included with the cabinet and must be screwed to the openings on the top section.

## 10. FIRE - DISPOSAL

### 10.1. OPENING THE CABINET AFTER THE FIRE

**CAUTION:** Depending upon the duration of the fire an ignitable vapour/air mixture may have formed; therefore remove all sources of ignition within a 10 metre radius of the cabinet before opening it.

- use only spark-free tools
- open the cabinets with extreme caution

### 10.2. DISPOSAL

The models can be disposed of once they have been dismantled and the materials sorted.

## 11. SAFETY CHECKS

### 11.1. IN GENERAL

As safety equipment the cabinets have to be checked for safety at least once per year. The next checking date can be taken from the check sticker on the outside of the door. This annual check can be carried out with the necessary care, and for securing your warranty claims in the case of fire, only by an authorised asecos employee (refer also to our service brochure regarding this).

**In addition, we recommend that you carry out a daily and monthly functional check:**

#### **Daily functional check**

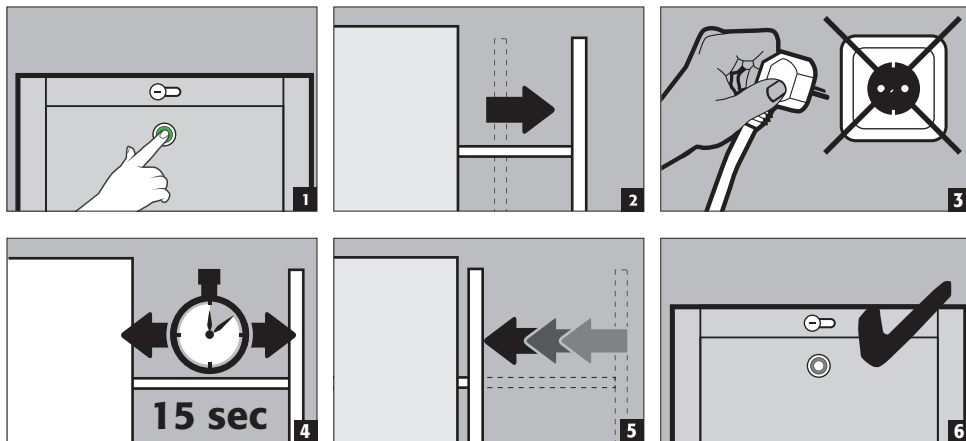
- applies to the bottom collecting sumps (collect and remove any leakages)

#### **Monthly functional check**

- perfect function of the drawers
- correct seating and condition of the fire protection seals

## 11.2. BATTERY CHECK

- Every six months after the maintenance check the battery should be checked by the user of the cabinet.

In case of damage please contact your dealer in order to have the cabinet repaired using original spare parts. The cabinets can be cleaned with a mild household cleaner and a soft cloth.

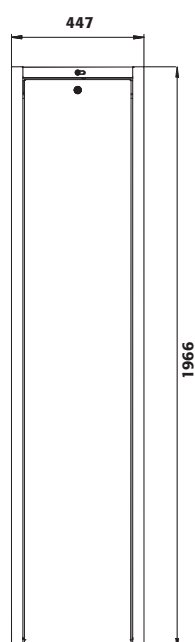
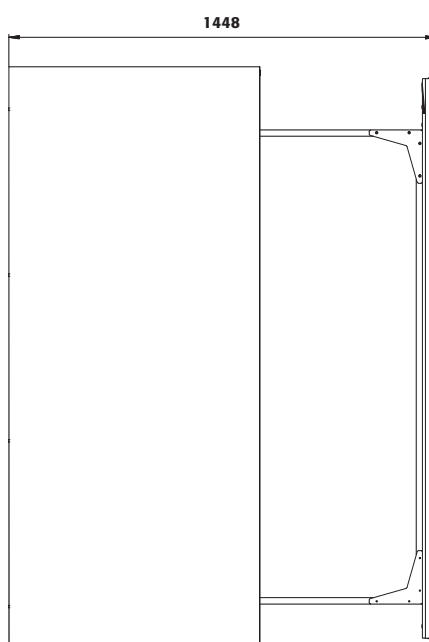
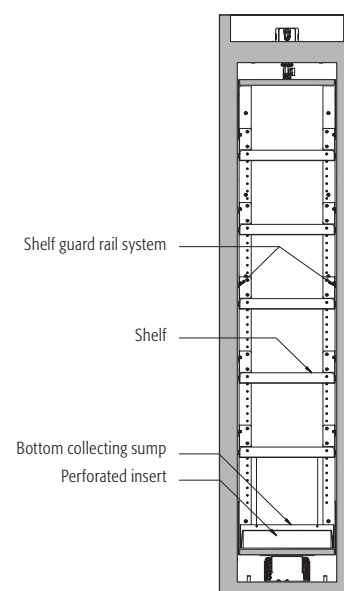
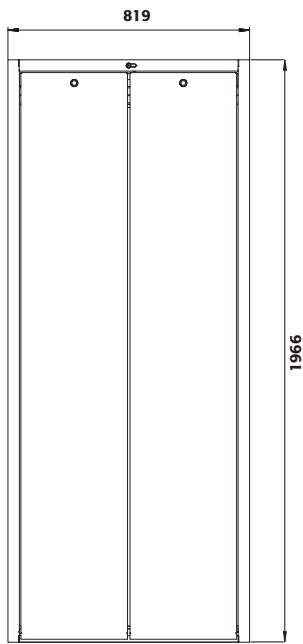
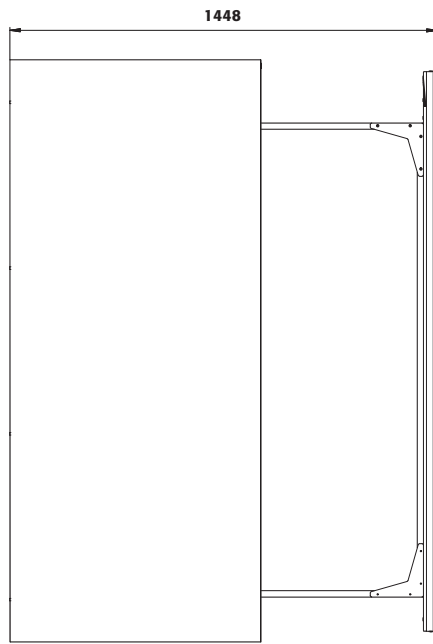
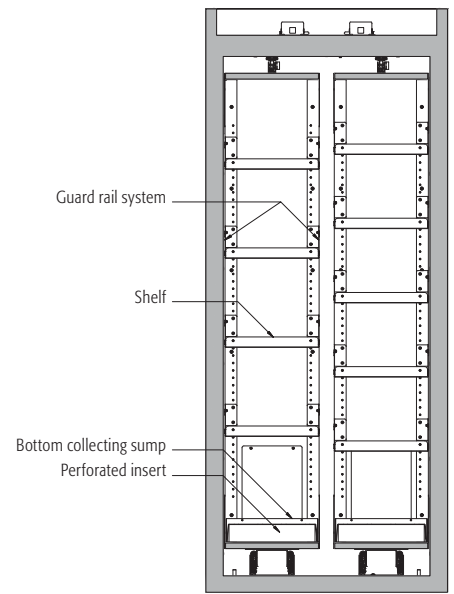
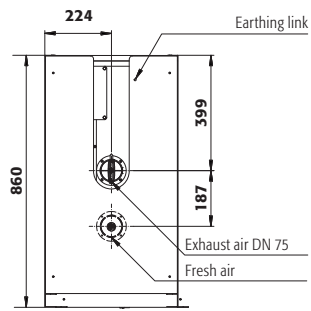
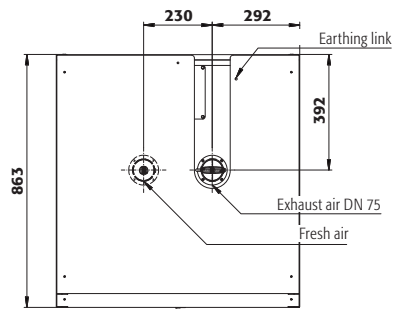
**CONTACT:** In the case of defects or complaints about our products (within and also after the warranty period), and for requesting safety checks or taking out a service contract, please contact our service hotline on: Tel: +44 1785 22 70-90 [info@asecos.co.uk](mailto:info@asecos.co.uk)

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