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User manual

**Extraction units**

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**Ventilatie-opzetunits**

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**Boîtiers de ventilation**

Manual de instrucciones

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---

**HFEA.8677 / HFEA.11442 / HFEA.17447 / HFEA.11570 / HFEA.15575 / HFEA.17448 / HFEA.8473 / HFEA.23866 / HFEA.24139**

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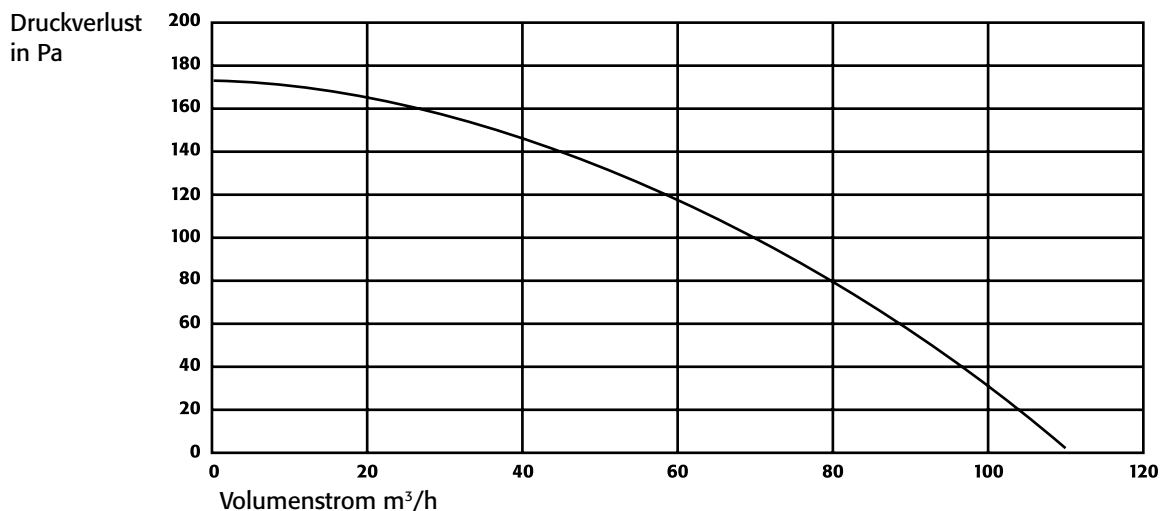
The extraction air unit is used for the technical extraction of air from the inside of cabinets. The extracted air stream does not get in contact with the electric current leading or metallic blank construction parts. The unit is suitable to extract cabinets which are defined in the EN 14470-1/2.

Extraction air unit: RL 94/9/EG (ATEX) und RL 73/23/EWG

 II 3/- G c IIC T4

## 2. Assembly

---

**HFEA.8677 / HFEA.11442 / HFEA.17447**

Install the extraction unit with the extraction air inlet on top of the cabinet. Connect the customer provided extraction air duct appropriate to the extraction air outlet of the unit. The extracted air from in side safety cabinet can either flow into a suitable ventilation system or must flow in safe place into the free.

**ATTENTION: Use plastic duct with cabinet models VBFA/VBFT/VBF type series 196.**

**HFEA.11570 / HFEA.15575 / HFEA.17448 / HFEA.23866 / HFEA.24139**

Ensure that all parts listed in the enclosed packing list are present.

**Wall consoles:** Install both triangular steel sheet parts with three screws each on the side of the extraction air ventilator housing.

**Choice of installation location:** Look for a location above or near the hazardous material cabinet, as close as possible to the outlet connector of the cabinet.

**Wall mounting:** Drill the necessary holes in the wall, remove the bore dust and insert the plastic rawlplugs. Attach the ventilator cowl, with the enclosed screws on the wall.

**HFEA.8473**

1. Bolt the enclosed brackets onto the side of the ventilation module.
2. Place the ventilation module with its inlet opening over the outlet connector of the underbench cabinet, and align it horizontally.
3. Mark the positions of the holes required for the brackets on the rear wall of the underbench cabinet. Spot-drill the marked holes with a small wood drill. Then position the ventilation module with the brackets in the appropriate positions, and fasten the brackets to the rear wall with the wood screws included.
4. Connect the outlet connectors (Ø 75 mm) on the ventilation module to the customer's exhaust air ducts. Note that the connection must exhibit a degree of elasticity.
5. Insert the mains plug into a socket that is earthed in accordance with regulations.

## 3. Put into operation

---

Put the power supply plug into a correctly grounded power socket.  
Activate the rocker switch (Position ON).

## 4. Extraction air monitoring

---

**HFEA.11442 / HFEA.15575 / HFEA.17447 / HFEA.17448 / HFEA.23866 / HFEA.24139**

Monitoring with integrated pressure leak detector. The green light turns off in case of failure.

## 5. Connections/repairs

---

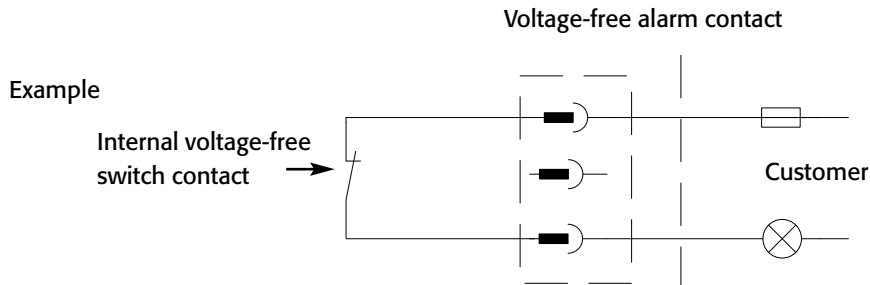
Ventilation connections as well as repairs shall only be done by specialists. In the event of damage the equipment has to be repaired or exchanged by the manufacturer.

## 6. Voltage-free alarm contact

### HF.EA.17447 / HF.EA.17448 / HF.EA.24139

The supplied three-pin plug is needed for using the voltage-free alarm contact. Only the two external terminals are connected to the plug! Internally a contact is switched (closed) as soon as the circulating air filter attachment reaches the prescribed volumetric flow.

Thus signalling is on hand in case of extraction air trouble or when your circulating air filter attachment is disconnected from the mains.

**The voltage-free alarm contact is not to be used for operating large loads. It is only envisaged for signalling a malfunction by means of a signalling lamp, horn and the like!**

## 7. Technical data

The figures in the table are provided as a guide, and are not to be used in the sense of assured properties.

### HF.EA.8677 / HF.EA.11442 / HF.EA.11570 / HF.EA.15575 / HF.EA.8473 / HF.EA.17447 / HF.EA.17448 / HF.EA.23866 / HF.EA.24139

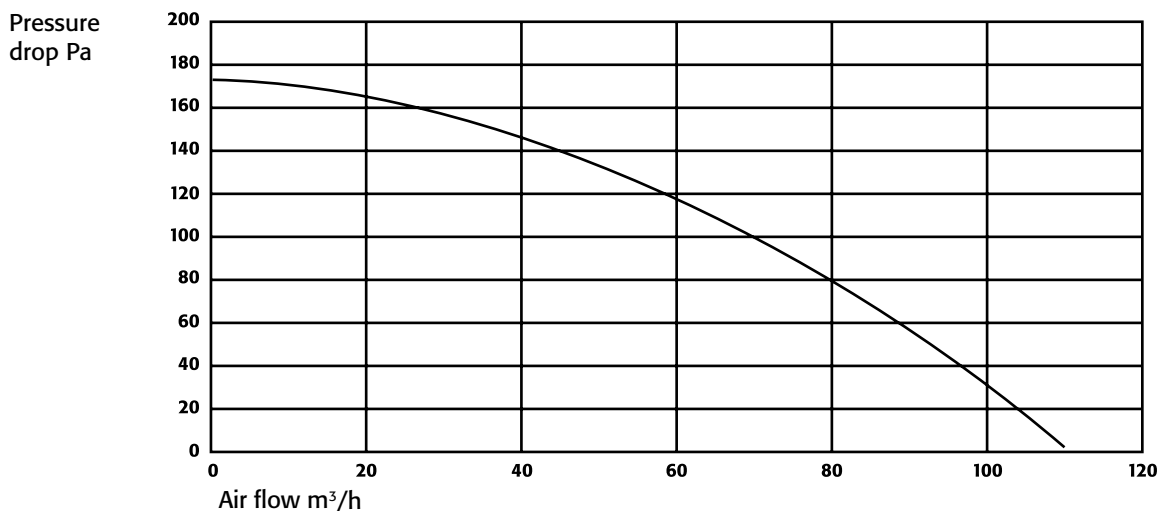
Dimensions W x D x H (mm)	200 x 400 x 200
Nominal rating	35 Watt
Nominal voltage	230 V
Frequency	50 Hz
Power consumption	0,15 A
Noise level	3m/35 dB (A)
Revolution Rpm	2450
Weight approx.	ca. 7 kg
Max. Length of extraction air duct by customer (including max. 2 pipe elbows)	15 m

### Load capacity of the voltage-free switch contact

Maximum switching-on current AC	230 V ~
Maximum switching-on current DC	48 V
Maximum permanent current	1A

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