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Umluftfilteraufsatz UFA.20.30

USER MANUAL

Circulating Air Filter UFA.20.30

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Recirculatie-opzetunit UFA.20.30

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TABLE OF CONTENTS

1. NOTES • GUIDELINES • GUARANTEE	8
1.1. General Safety Notes	8
1.2. Guarantee	8
1.3. Intended Application	8
1.4. Development and Manufacture	9
2. DESIGN	9
2.1. Material	9
2.2. Accessories	9
3. MODE OF OPERATION	9
3.1. Function	9
3.2. Display	10
4. START UP AND OPERATION	10
4.1. Start up	10
4.2. Self-test	10
5. MALFUNCTIONS • WARNINGS	11
6. VENTILATION ADJUSTMENT • MONITORING	11
7. VOLTAGE-FREE ALARM CONTACT	11
8. TECHNICAL DATA	12
9. REPAIRS	12
10. FUNCTIONAL CHECK • CARE • MAINTENANCE	12
11. VIEWS • CROSS-SECTION	13
12. ATEX	13

1. NOTES • GUIDELINES • GUARANTEE

1.1. GENERAL SAFETY NOTES

- Take notice of the legislation and provisions to be applied in dealing with hazardous materials.
- Only use the extraction air inlet on the top in a proper condition.
- Take note of all information in these operating instructions.
- Follow the instructions of the safety specialists, technical monitoring services (BG, GAA, pharmaceuticals councils), construction supervision etc.
- Observe accident prevention regulations and workplaces ordinances.
- Regularly examine the extraction air inlet on the top as to proper functioning (check on the LED displays for extraction air volumetric flow and filter saturation).

1.2. GUARANTEE

The guarantee for this product is agreed between you (the customer) and your dealer (the seller). Independently of this, the manufacturer provides a guarantee for the extraction air inlet on the top for 24 months after the date of sale. This device is classified as safety-relevant equipment and must be checked at least every year, or at shorter intervals depending on load. Inspections and filter changes must be carried out by specialised staff authorised by the manufacturer, using original spare parts from the manufacturer. If this inspection is not carried, or not carried out by specialised staff authorised by the manufacturer, the customer's guarantee claims against the manufacturer will be invalidated.

1.3. INTENDED APPLICATION

The purpose of the appliance is to ventilate from ex-zone 2 or the non ex-zone of cabinets. The extraction air inlet on the top is especially quiet and thus ideal for using in the immediate working environment. Suitable for cabinet models **S-CLASSIC-90 / S-PEGASUS-90** (formerly VBF/T/A.196.120/90/60/6.7 and VBF.129.120/60), **Q-LINE** (formerly VBF.195.120/.90/.60) together with attachment adapter HFA.15776: **XL-LINE** drum cabinets (formerly VBF.221.108/.153) and **S-CLASSIC-30** (formerly VBF.197.xxx).

**CAUTION:**

Installation may only take place in a non-ex-zone!

The extraction air is cleaned to 99.999 % of hydrocarbons by the integrated activated carbon filter so that it can be directly given off to the ambient air. The extraction air flow containing pollutants does not come into contact with energised components. Hence the appliance is suitable for suctioning out cabinets defined in EN 14470 Part 1 resp. DIN 12925 Part 1. There is thus no need for an extraction air line leading to a safe place outdoors.

A CE conformity attestation (also in acc. with EC directive 94/9/EC - ATEX directive) is on hand.

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
SN: The type plate with the serial number is situated on the outside of the housing of the UFA.20.30.

EN

2. DESIGN

2.1. MATERIAL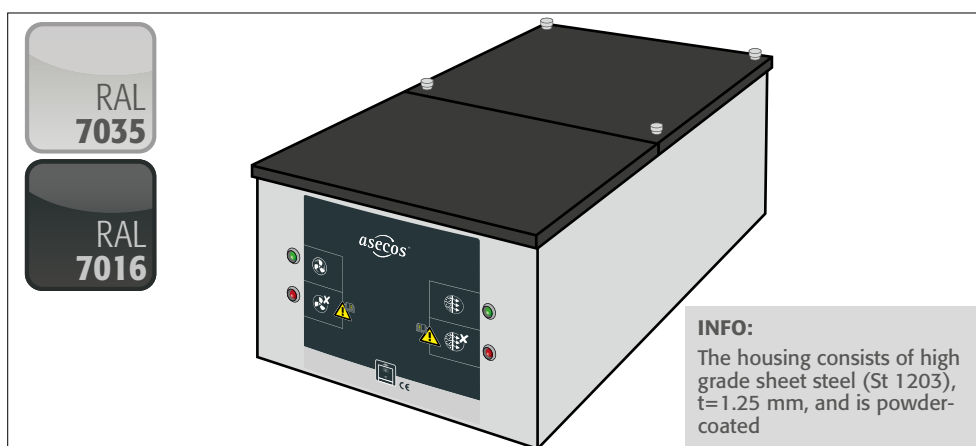


2.2. ACCESSORIES

Activated carbon filter (3,500 g) incl. suspended solids preliminary filter.

3. MODE OF OPERATION

3.1. FUNCTION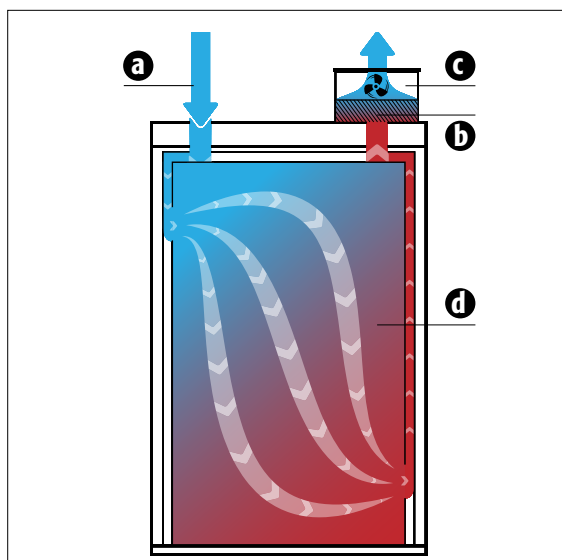

- a** fresh air
- b** activated carbon filter
- c** ventilator
- d** gases and vapours in cabinet

CAUTION:


Filter can only retain solvent vapours (hydrocarbons). Please consult us if other materials are to be stored.

Safe and faultless function is shown by the ventilation and filter monitoring mechanism which is fitted as standard. (3.2)

3.2. DISPLAY

MALFUNCTION:
for actions see
6. Malfunctions




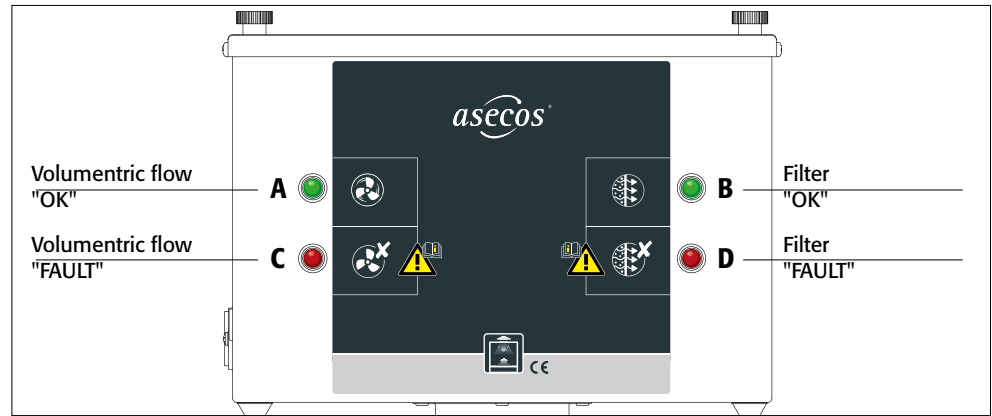


4. START UP AND OPERATION

4.1. START UP







CAUTION:

Observe the notes in item
1.3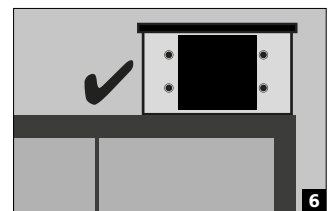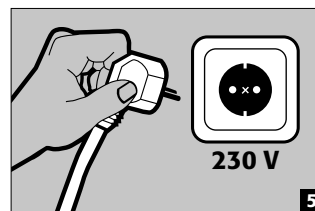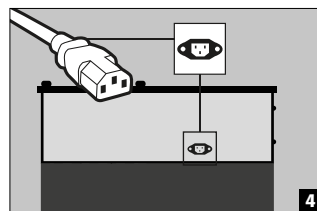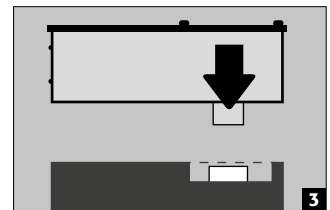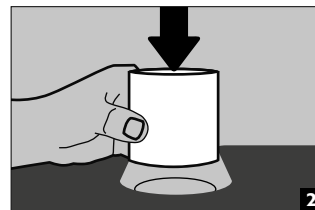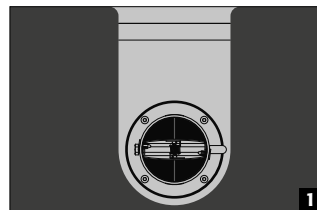



4.2. SELF-TEST

After inserting the mains plug, the device performs a self-test: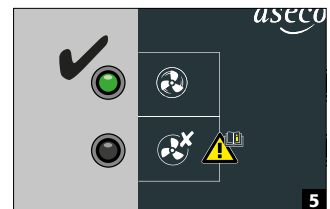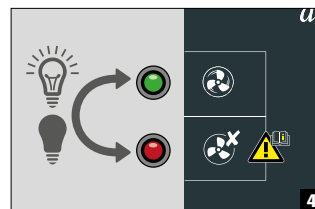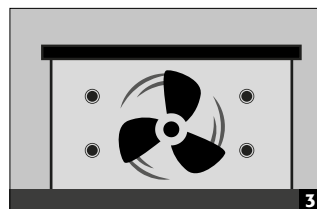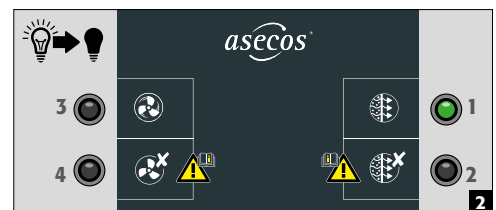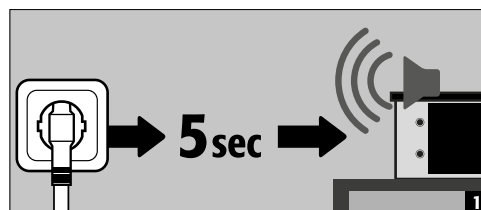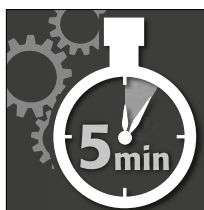

CAUTION:

The device is designed for
continuous operation!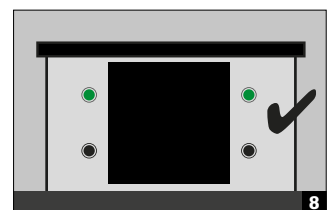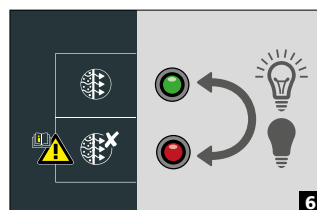

5. MALFUNCTIONS - WARNINGS

FAULT	REASON – ACTION	NOTES - must be observed!
1 Inserting the mains plug <ul style="list-style-type: none"> • A and B are not illuminated • Ventilator is off 	1. No current in the mains socket or 2. effective fuse in fullymains plug is inserted insulated socket or 3. defective energy saving lamp illuminant.	The C13 cable must be removed before replacing the fuses
2 after the self-test <ul style="list-style-type: none"> • C and D flash • 5 beep sounds • Ventilator is off 	1. Possible defect of the device or 2. Defect of the monitoring circuitry	please contact your dealer or our costumer service: +49 6051 92 20 0* (* at local rates in the german fixed network)
3 during operation <ul style="list-style-type: none"> • A and C are illuminated • intervall sound (30 sec.) • Ventilator is off 	1. Foreign matter may possibly be in or on the (air flow) extraction air opening (to be removed) or 2. Possible defect of the integrated extraction air ventilator.	please contact your dealer or our costumer service: +49 6051 92 20 0* (* at local rates in the german fixed network)
4 during operation <ul style="list-style-type: none"> • A is illuminated • D flashes 	The filter has almost reached its saturation limit; a timely (filter monitoring) replacement is called for!	please contact your dealer or our costumer service: +49 6051 92 20 0* (* at local rates in the german fixed network)
5 during operation <ul style="list-style-type: none"> • A and D are illuminated • Alarm tone 2 sounds out every hour • Ventilator is off 	The filter is almost fully saturated and the ventilator has been switched off. The filter must now be changed quickly.	please contact your dealer or our costumer service: +49 6051 92 20 0* (* at local rates in the german fixed network)
6 during operation <ul style="list-style-type: none"> • A and D are illuminated • Alarm tone 1 sounds out every hour 	The permissible operating time of the filter has expired. The filter must now be changed quickly.	please contact your dealer or our costumer service: +49 6051 92 20 0* (* at local rates in the german fixed network)

ALARM SOUNDS

There are two different alarm sounds. If both are simultaneously activated, they ring out after one another at an interval of approx. 6 seconds.

Tone sequence 1: short – short – long

Tone sequence 2: long – short – long – short

6. VENTILATION ADJUSTMENT - MONITORING

The ventilator capacity is sufficient for a minimum ten-fold change of air for those cabinet models listed under the „Purpose“ item.


The faultless ventilation function is automatically controlled by a monitoring electronics system.

A differential pressure measurement device monitors air volumes in the extraction air shaft.

An alarm is triggered (**C3**) when the air rated value is undercut by around 50 %. The optical alarm only stops when the rated value is again attained.

7. VOLTAGE-FREE ALARM CONTACT

The supplied three-pin plug is needed for using the voltage-free alarm contact. Only the two external terminals are connected to the plug! Internally a contact is switched (opened) as soon as the extraction air inlet on the top reaches the prescribed volumetric flow. An extraction air fault or filter saturation can therefore be signalled in this way. Furthermore it is signalled when the extraction air inlet on the top is disconnected from the mains.

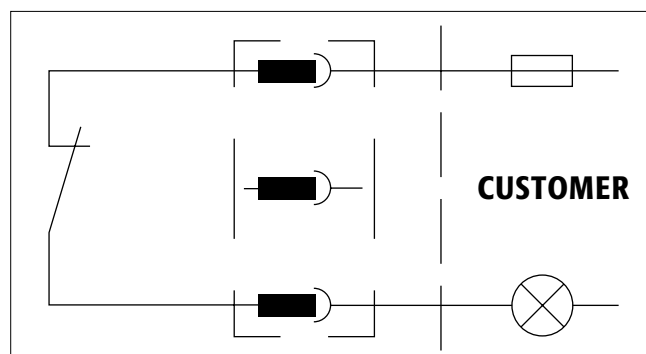CAUTION:

The voltage-free alarm contact is not to be used for operating large loads.

It is only envisaged for signalling a malfunction of the extraction air inlet on the top by means of a signalling lamp, horn and the like.

8. TECHNICAL DATA

Technical data	UFA.20.30
External dimensions (H×W×D mm)	210×300×550
Nominal power	30 Watt
Nominal voltage	230 V ~± 10%
Frequency	50 Hz
Current consumption	0.13 A
Noise level	39 dB
Net weight (without Filter)	12 kg
Activated carbon filter weight (inkl. Füllung 3500 g)	4.2 kg
Volumetric flow (mounted on the cabinet)	25 m³/h
Load capacity of the voltage-free alarm contact	
Max. nominal voltage AC	250 V ~
Max. permanent current at 250 V AC	6 A
Max switching power (ohms)	1500 VA
Max switching power (elektro-magnetic)	300 VA
Single-phase motor load (inductive)	185 Watt
Max. switching current at 30 V DC	6 A
Max. switching current at 115 V DC	0.2 A
Max. switching current at 230 V DC	0.12 A

CAUTION:

The figures in the table are provided as a guide, and are not to be used in the sense of assured properties. They conform to the state of the art as and when the operating instructions were printed. Changes to these values without notice as a result of further technical development are explicitly permitted.

9. REPAIRS

The extraction air inlet on the top is only to be repaired by specialists trained for this very purpose. Given damage the appliance is to be repaired or replaced by the manufacturer. Please contact the customer service.

10. FUNCTIONAL CHECK - CARE - MAINTENANCE

To ensure reliability, the extraction air inlet on the top should be serviced and checked as to functioning by a specialist at least once a year. At the same time the fitted filter unit is to be replaced at least once a year - or sooner should the saturation limit be reached.

The integrated ventilation and filter saturation monitoring system electronically checks on functionality. As a result, the extraction air inlet on the top function is constantly examined during operation. The activated carbon filter of the extraction air inlet on the top must be replaced soon after the red LED of the filter display blinks. Only our trained service technicians are to change the filter, which is to be rapidly requested under the Service hotline. Having the filter changed by the asecos service ensures that contaminated activated carbon is recycled in a defined disposal chain. Clean the outside of the extraction air inlet on the top with a soft cloth and a common, mild household cleaner. Before cleaning, disconnect the extraction air inlet on the top from the mains by pulling out the C13 cable! Make sure that no moisture can penetrate into the inside of the extraction air inlet on the top during cleaning.


SERVICE: In case of questions please contact Customer Service!

Tel: +49 6051 92 20 0* (* at local rates in the german fixed network)

11. VIEWS - CROSS-SECTION

- Ventilator **1**
 Gas sensor **2**
 Filter **3**
 Voltage-free alarm contact **4**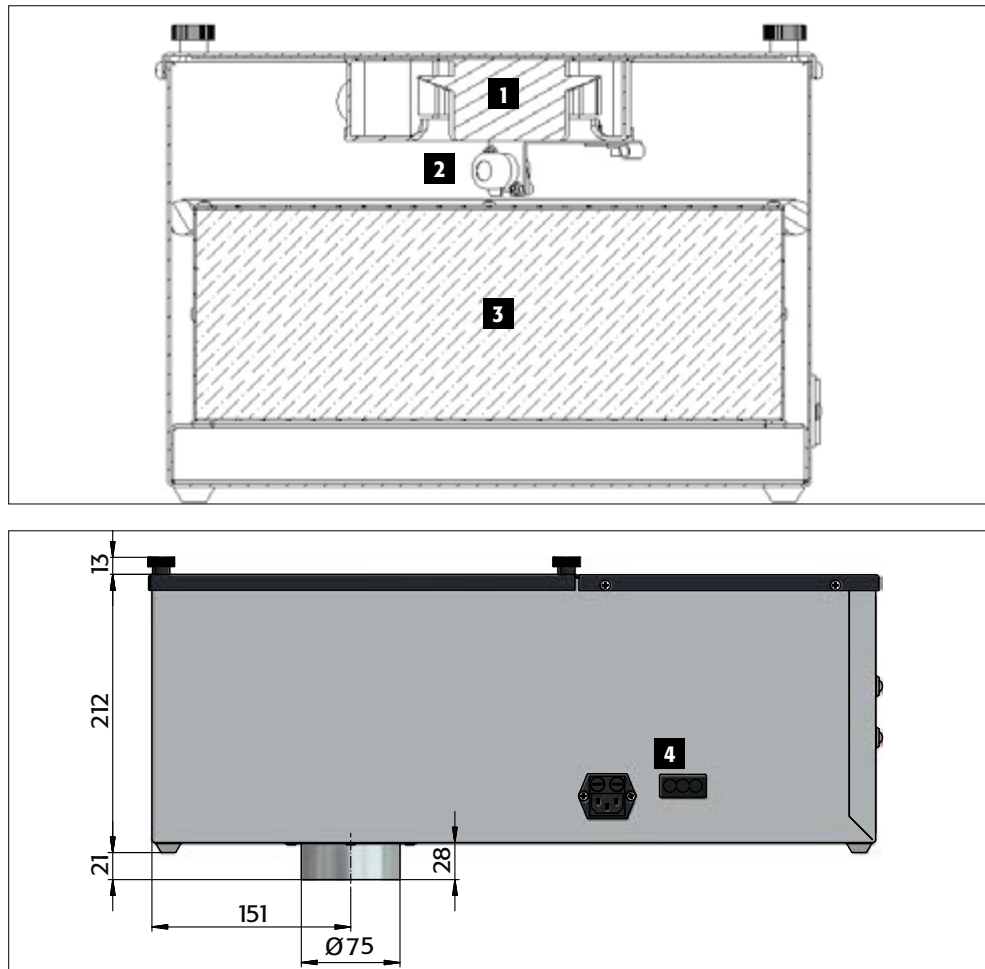

12. ATEX

	II	3/- G	Ex ic nA	IIB	T4	Gc
1	2	3	4	5	6	7

1	Marking of explosion-proof devices
2	Device group in accordance with directive 94/9/EC (Atex 95) II non-underground operation
3	Device categories of the assemblies used 3/- extraction from explosion zone 2, installation exclusively in non-explosive zones G Gas
4	Ignition protection classes Ex ic intrinsic safety with the protection level "ic" – not capable of causing ignition in undisturbed operation nA non-sparking
5	Explosion groups in accordance with EN 60079-0:2011 IIB the group has a safe gap value (MESG) of 0.5-0.9 mm, suitable e.g. for use of ethylene
6	Temperature class T4 max. surface temperature of the equipment 135 °C
7	Protection level EPL pursuant to IEC 60079-0 (2007) 3/- G Gc Suitability for extraction from explosion zone 2

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asecos GmbH